

AppliChrom® ABOA SugarSep-Oligo



28/09/2015 Application Note

Oligosaccharidanalytics in Water



Total chromatogram

AppliChrom ABOA SugarSep-Oligo, 300x8mm
"Maltodextrin 19", Eluent: H₂O, 0,25ml/min, 70°C



Detail enlargement

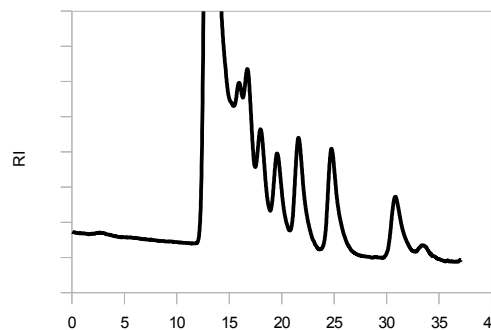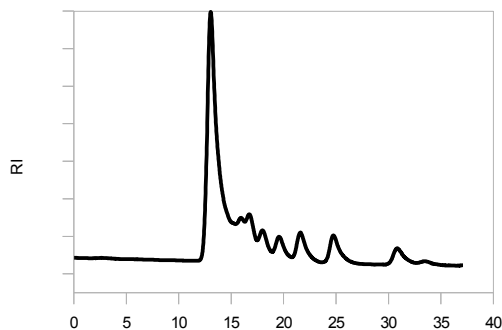

AppliChrom ABOA SugarSep-Oligo, 300x8mm
"Maltodextrin 12", Eluent: H₂O, 0,25ml/min, 70°C

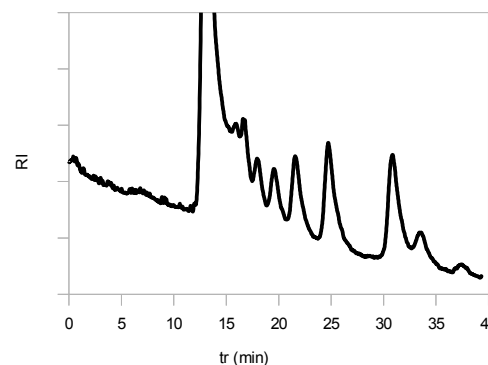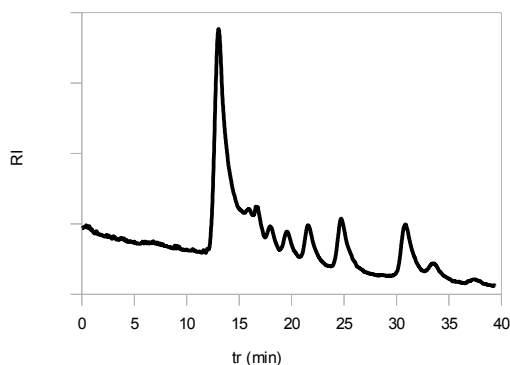

AppliChrom ABOA SugarSep-Oligo, 300x8mm
"Maltodextrin 6", Eluent: H₂O, 0,25ml/min, 70°C

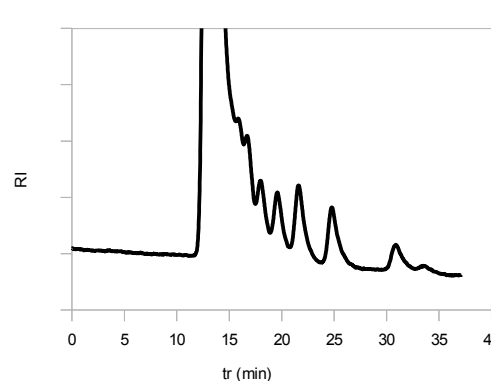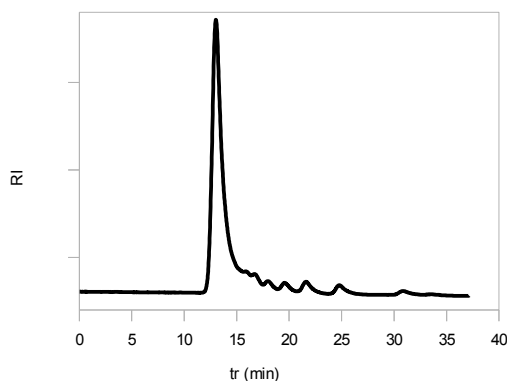

Order Information:

SASOL101508	HPLC-Column AppliChrom® ABOA SugarSep-Oligo, 10µ 150mm x 8mm	Price available upon request
SASOL103008	HPLC-Column AppliChrom® ABOA SugarSep-Oligo, 10µ 300mm x 8mm	Price available upon request
SASOL10508	Precolumn AppliChrom® ABOA SugarSep-Oligo, 10µ, 50mm x 8mm	Price available upon request

AApplication note 09/2015, changes and errors reserved. Not intended for pharmaceutical, diagnostic or medical use, for research and laboratory only. Only terms and conditions of AppliChrom are valid.