

Our new electronic crimpers and decappers for 8mm, 11mm, 13mm and 20mm crimp caps fit standard laboratory vials and seals with aluminum or two-part aluminum/steel caps. The electronically adjustable cycle is accurate and reproducible and can be adjusted with the control pad on top of the unit. The tools are ergonomically designed to reduce strain and arm injury associated with many manual crimping tools.

- Powerful motor with fast cycle
- Longer lasting battery with Lithium Ion technology
- Advanced charging system including warning light when recharging is necessary
- Can be operated while charging
- Convenient controls on top of unit



Quick Start

1. Step - Plug in and charge until LED shows permanent green light




2. Step - Perform a crimp test before using your sample vials



Adjust Crimper

Too Tight

Too Loose

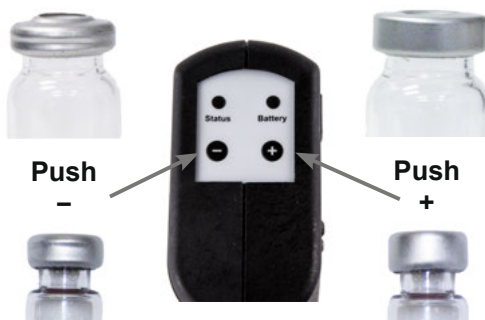

Adjust Decapper

Not Enough



Note: Use the Electronic Crimper for Aluminium Crimp Caps only. For Magnetic Caps use the HPS Crimp Tool.