

EVOKE™

C18 Columns

High performance reversed-phase separations at any scale


- High efficiency columns designed to resolve complex samples
- Unique gas-phase endcapping for added hydrophobicity and inertness
- Seamless scalability with particles from 1.8 μm to $>20 \mu\text{m}$
- Proprietary packing technology ensures column to column reproducibility




Scalable Separations across a Wide Range of Particle Sizes

Separation of NSAIDs on Evoke C18



100 x 2.1 mm; 1.8 μm
Flow Rate: 0.4 mL/min

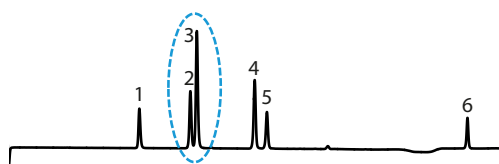

100 x 4.6 mm; 3 μm
Flow Rate: 2.0 mL/min

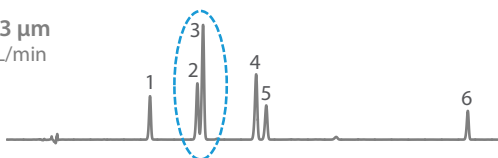

100 x 4.6 mm; 5 μm
Flow Rate: 2.0 mL/min

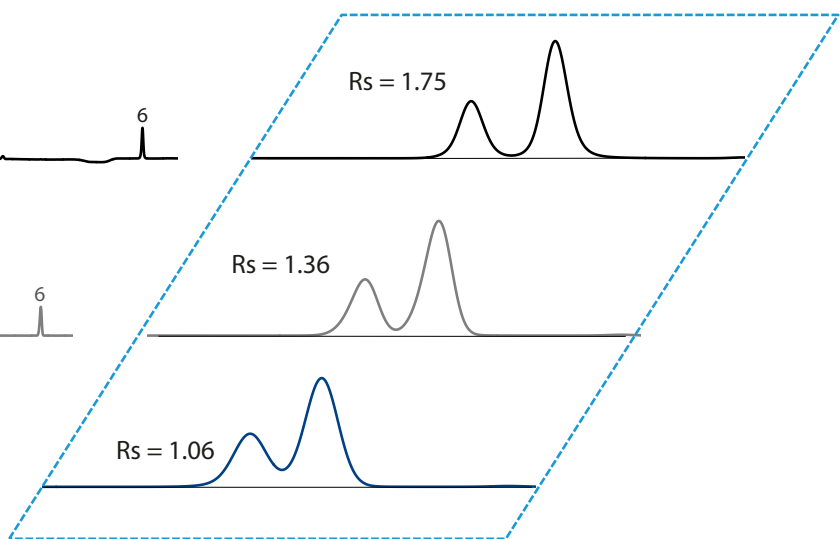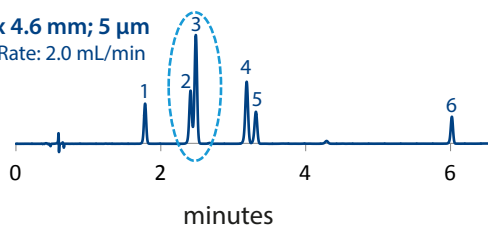

Peak IDs: 1) Indoprofen 2) Ketoprofen 3) Naproxen 4) Fenoprofen 5) Flurbiprofen 6) Ibuprofen

Mobile Phase: A) H_2O + 0.1% TFA; B) ACN + 0.1% TFA

System: Shimadzu Nexera UC

Detection: UV 220 nm


Gradient:

Time (min)	%B
0	40
4	70
6	95

A new line of C18 columns, optimized for high efficiency separations and excellent column to column reproducibility, improves productivity and solves complex separation challenges. Evoke™ C18 columns are packed with high purity silica and endcapped using a unique gas-phase technology. Gas-phase endcapping leads to improved surface coverage, shielding acidic silanol groups for a highly inert, hydrophobic surface that results in excellent separation kinetics and provides unique selectivity in comparison to other C18 phases. Additionally, it allows the use of higher concentrations of organic mobile phase for greater sensitivity by electrospray ionization mass spectrometry and improved loading capacity for preparative applications.

High efficiency separations can be achieved on fully porous particles, ranging from 1.8 μm (analytical scale) to greater than 20 μm (process scale purifications). Our most popular column sizes are stocked for immediate delivery.

High Capacity C18 Phase

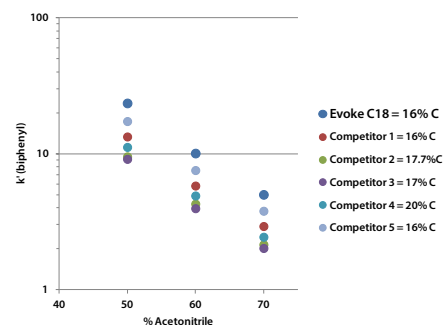

Unique gas-phase endcapping promotes hydrophobic retention for unique selectivity and greater retention over leading high carbon load phases

EVOKE™ C18 Columns

ORDERING INFORMATION

1.8 µm

DIMENSIONS	CATALOG #
5 cm x 2.1 mm	1-760106-300
10 cm x 2.1 mm	1-760107-300
2 cm x 3 mm	1-760103-300
5 cm x 3 mm	1-760104-300
10 cm x 3 mm	1-760105-300
5 cm x 4.6 mm	1-760102-300
10 cm x 4.6 mm	1-760101-300

3 µm

DIMENSIONS	CATALOG #
5 cm x 2.1 mm	1-760206-300
10 cm x 2.1 mm	1-760207-300
15 cm x 2.1 mm	1-760209-300
2 cm x 3 mm	1-760203-300
5 cm x 3 mm	1-760204-300
10 cm x 3 mm	1-760205-300
15 cm x 3 mm	1-760210-300
5 cm x 4.6 mm	1-760202-300
10 cm x 4.6 mm	1-760201-300
15 cm x 4.6 mm	1-760208-300
25 cm x 4.6 mm	1-760211-300

5 µm

DIMENSIONS	CATALOG #
3 cm x 2.1 mm	1-760319-300
5 cm x 2.1 mm	1-760306-300
10 cm x 2.1 mm	1-760307-300
15 cm x 2.1 mm	1-760309-300
5 cm x 3 mm	1-760304-300
10 cm x 3 mm	1-760305-300
3 cm x 4.6 mm	1-760318-300
5 cm x 4.6 mm	1-760302-300
10 cm x 4.6 mm	1-760301-300
15 cm x 4.6 mm	1-760308-300
25 cm x 4.6 mm	1-760311-300
25 cm x 10 mm	1-760312-300
10 cm x 21.1 mm	1-760303-300
15 cm x 21.1 mm	1-760310-300
25 cm x 21.1 mm	1-760315-300
15 cm x 30 mm	1-760316-300
25 cm x 30 mm	1-760313-300
15 cm x 50 mm	1-760317-300
25 cm x 50 mm	1-760314-300

10 µm

DIMENSIONS	CATALOG #
5 cm x 2.1 mm	1-760406-300
5 cm x 4.6 mm	1-760402-300
10 cm x 4.6 mm	1-760401-300
15 cm x 4.6 mm	1-760407-300
25 cm x 4.6 mm	1-760411-300
25 cm x 10 mm	1-760412-300
15 cm x 21.1 mm	1-760410-300
25 cm x 21.1 mm	1-760409-300
50 cm x 21.1 mm	1-760404-300
25 cm x 30 mm	1-760413-300
50 cm x 30 mm	1-760405-300
25 cm x 50 mm	1-760403-300

Evoke™ C18 Media Specifications

Carbon Content:	16%
Pore Size:	100Å
Surface Area:	~320 m ² /g
pH Range:	2 - 8
Particle Sizes:	1.8, 3, 5, 10 µm and greater*

*Particles larger than 10 µm available upon request

Contact Regis today to discuss how our reversed-phase columns can improve your separations.

BGB gc|Lc
MS|CE

www.bgb-shop.com

Switzerland: BGB Analytik AG • Rohrmattstrasse 4 • 4461 Böckten • Phone +41 61 991 00 46 • Fax +41 61 991 00 25 • sales@bgb-analytik.com

BGB Analytik SA • Route de Pré-Bois 20 • 1215 Genève 15 • Phone +41 22 788 49 43 • Fax +41 22 788 49 45 • sales.fr@bgb-analytik.com

Benelux: BGB Analytik Benelux B.V. • Daltonstraat 17 • 3846 BX Harderwijk • Phone +31 341 700270 • Fax +31 341 700271 • sales.benelux@bgb-analytik.com

France: BGB Analytik France S.A.S. • 81 Vie de l'Etraz • 01630 St. Jean de Gonville • Phone +33 450 488567 • Fax +33 450 562378 • sales.fr@bgb-analytik.com

Germany: BGB Analytik Vertrieb GmbH • Josefstrasse 19a • 79618 Rheinfelden • Phone +49 7623 7173110 • Fax +49 7623 7173199 • sales.de@bgb-analytik.com